



CGQ / BioR Customized

ONLINE BIOMASS MONITORING FOR PHOTOTROPHIC APPLICATIONS

The CGQ and CGQ BioR allow for non-invasive online biomass monitoring in shake flasks and bioreactors.

Cell Growth Quantifier (CGQ) and CGQ BioR



A CGQ system consists of three components: the sensor(s), the hub, and the DOTS Software.

Hardware & Software Components

CGQ Sensor/ CGQ BioR sensor




The CGQ sensor plates are positioned under the shake flask, the BioR sensors are attached to the wall of the bioreactor. Both are designed to measure the biomass continuously and non-invasively.

Hub



The hub bundles the data from all monitored flasks or reactors and sends it to the DOTS Software.

DOTS Software

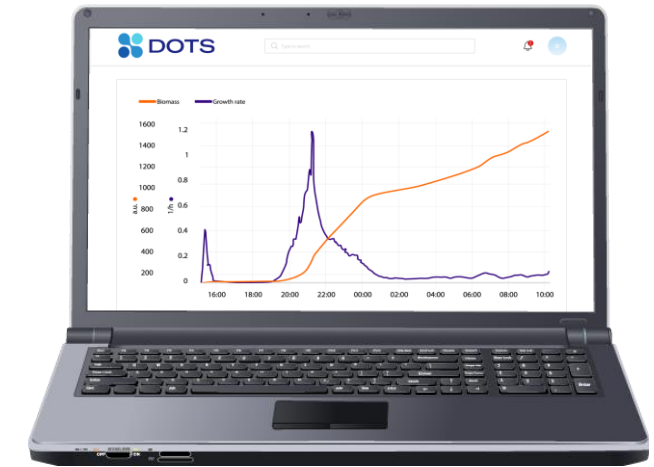

DOTS Software enables a simplified control of sensors and visualizes the received data from all monitored vessels in real-time.

The DOTS Software enables easy sensor handling and experiment set up.

Exemplary Screenshots

Create an experiment with pre-defined application templates

DOTS Type to search

1 Basic Settings 2 Device Assignment

Basic information

Quick start application templates Custom application templates

Biomass-based feeding (S... Select template

Enable template configuration step

Experiment name *
MET25 induction strain A

Number of objects *
1

Project *
Prosugar

Advanced settings

MET25 induction strain A

Biomass-based feeding (Shake flask)

Feeding (LIS) Planned
No device connected

Biomass monitoring (CGQ) Planned
No device connected

OD600 (Offline) Planned

Assign sensors to planned experiments via drag and drop

DOTS Type to search

Basic Settings General Task Configuration Replicate configuration Device Assignment

Filter

Type to search

Drag and drop device in order to connect it to process / task.

LIS-0025478
LIS-LIS0-254869 / 25
80% Not connected Unassigned

LIS-00-25478

CGQ-SP-02548
CGQ-8-0025 / Port1
Not connected Unassigned Free

MET25 induction strain A (0/1) (0/1)

Biomass-based feeding (Shake flask)

Feeding (LIS) Planned
No device connected

Biomass monitoring (CGQ) Planned
No device connected


OD600 (Offline) Planned

Exit wizard Back Create


The DOTS Software provides a comprehensive overview of your experiments and visualizes your data in real-time.

DOTS Software Modules for Data Visualization

The dashboard schema provides an overview over all running, planned, or finished experiments

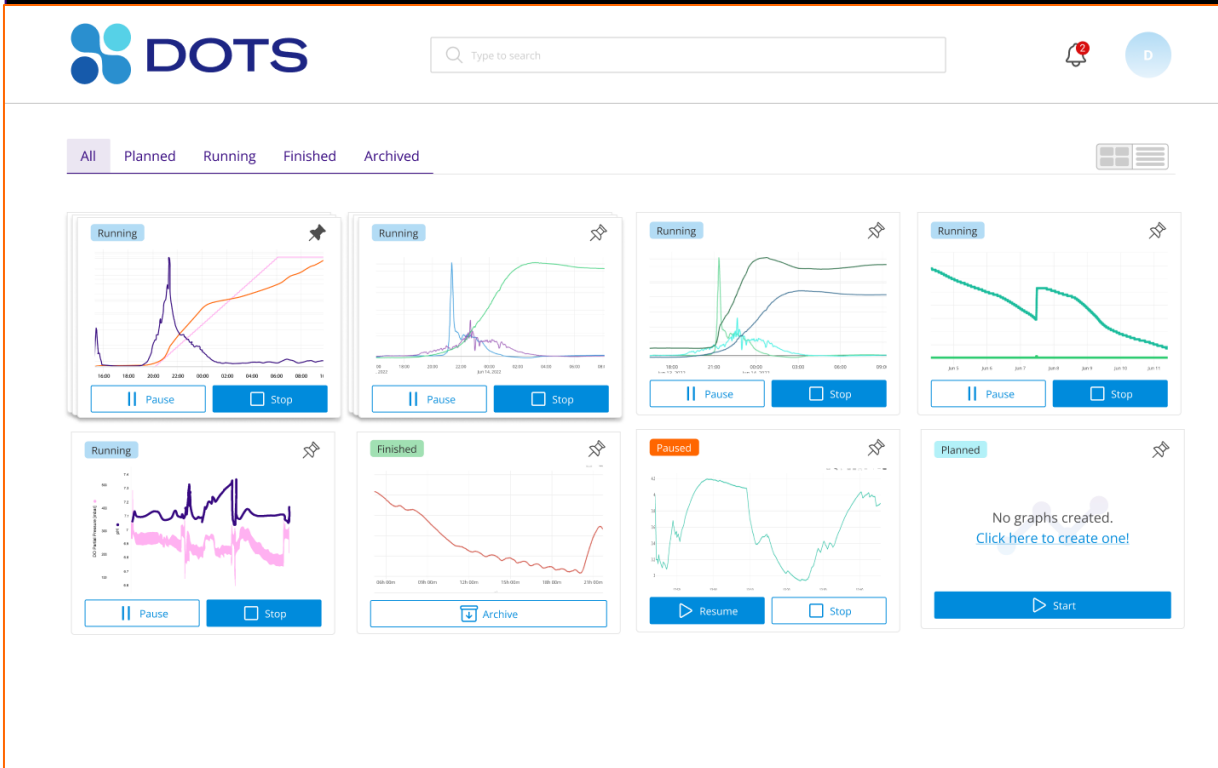


Data visualization tools enable a customized display of data in real-time

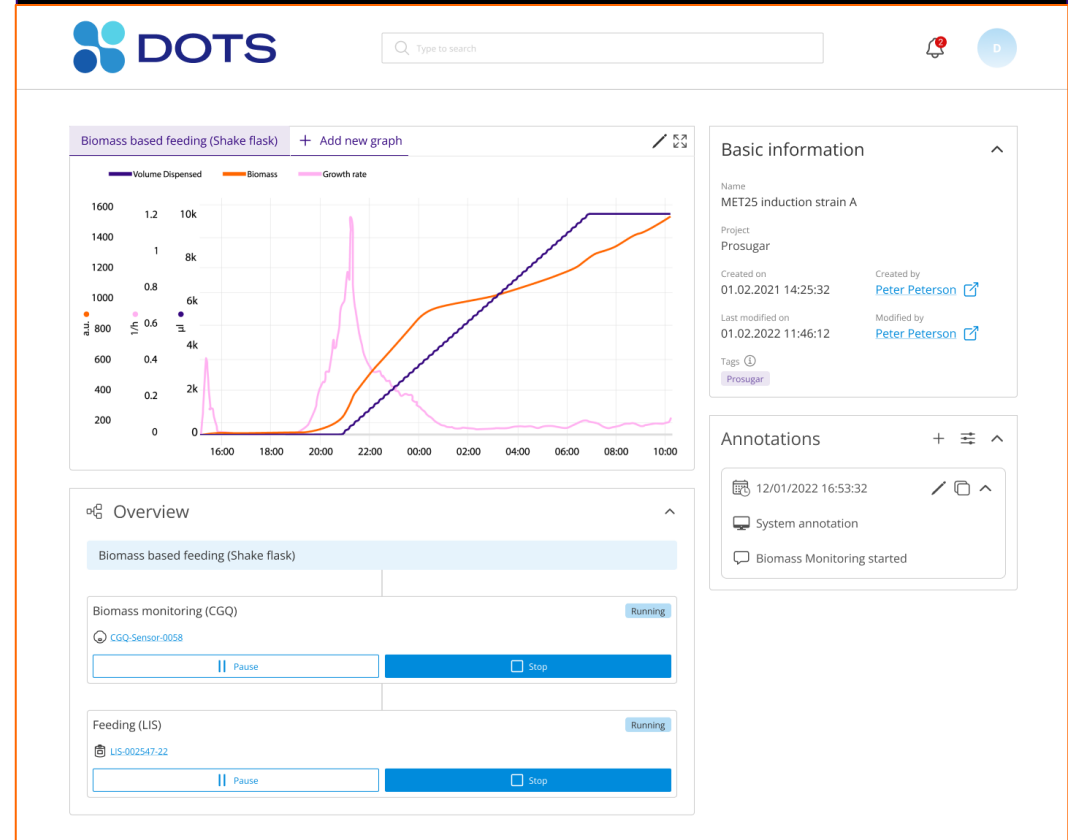


The CGQ can be installed in any phototrophic shaker and be used with any available flask size.

Exemplary Picture of CGQ Sensors Mounted in a Phototrophic Shaker

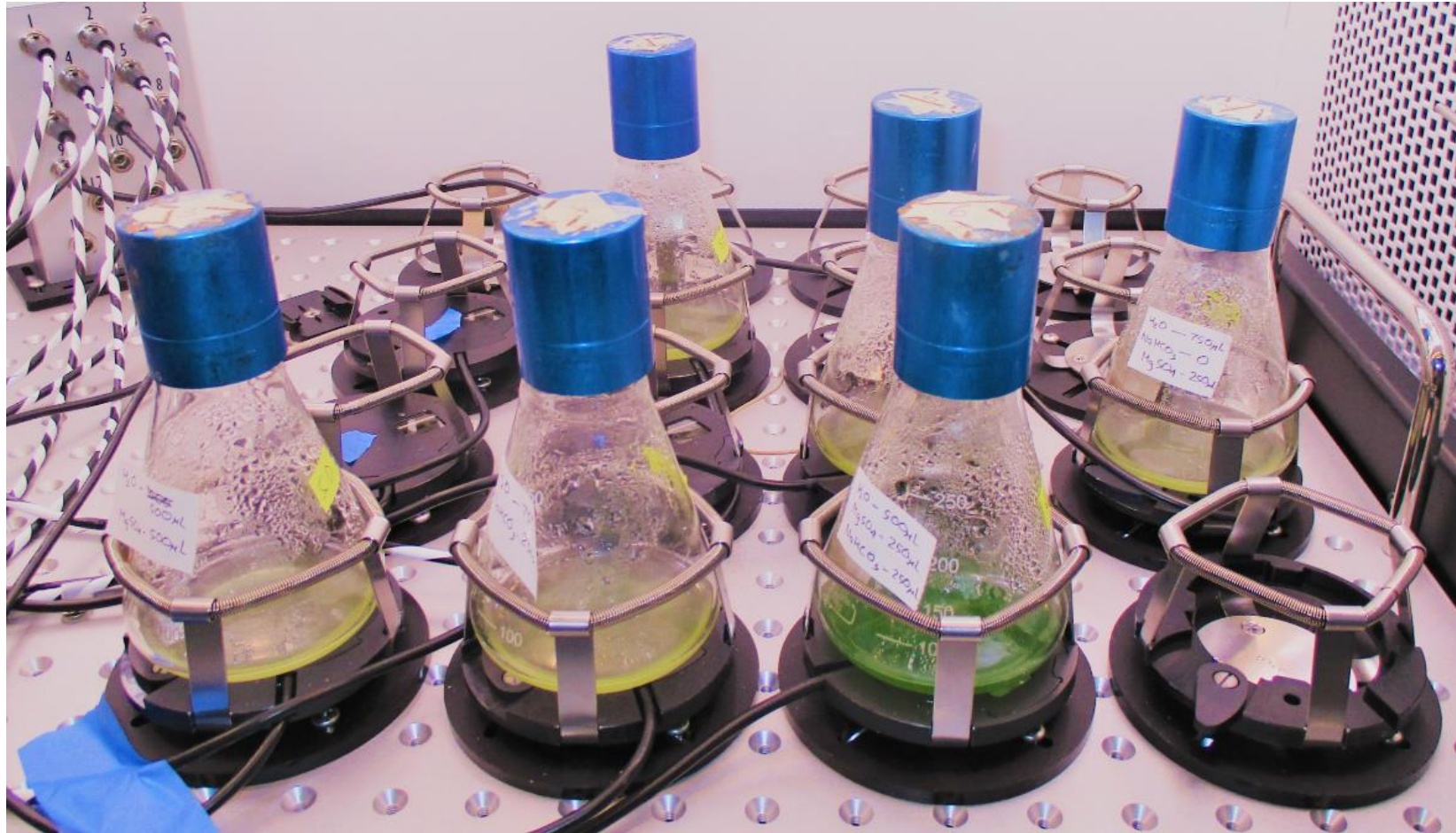


The CGQ BioR is perfectly suitable for specialized phototrophic bioreactors (photobioreactors) such as tubes or bags.

Exemplary Pictures of CGQ BioR Sensors Mounted to Different Types of Phototrophic Reactors

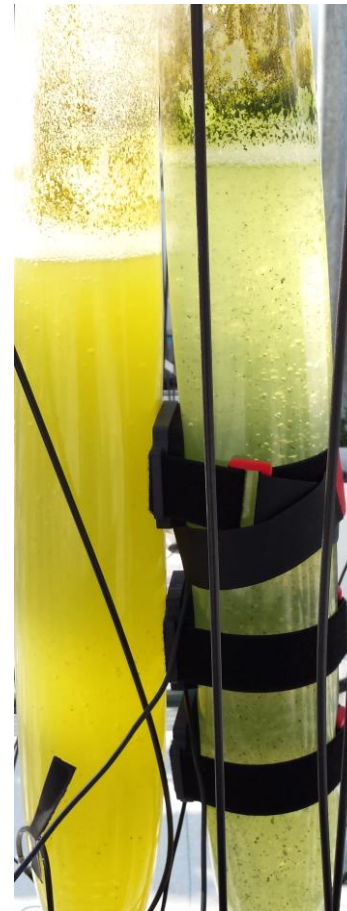


CGQ and CGQ BioR use backscattered light measurements allowing for accurate biomass monitoring in shaken cultures and bioreactors.

Measurement Principle

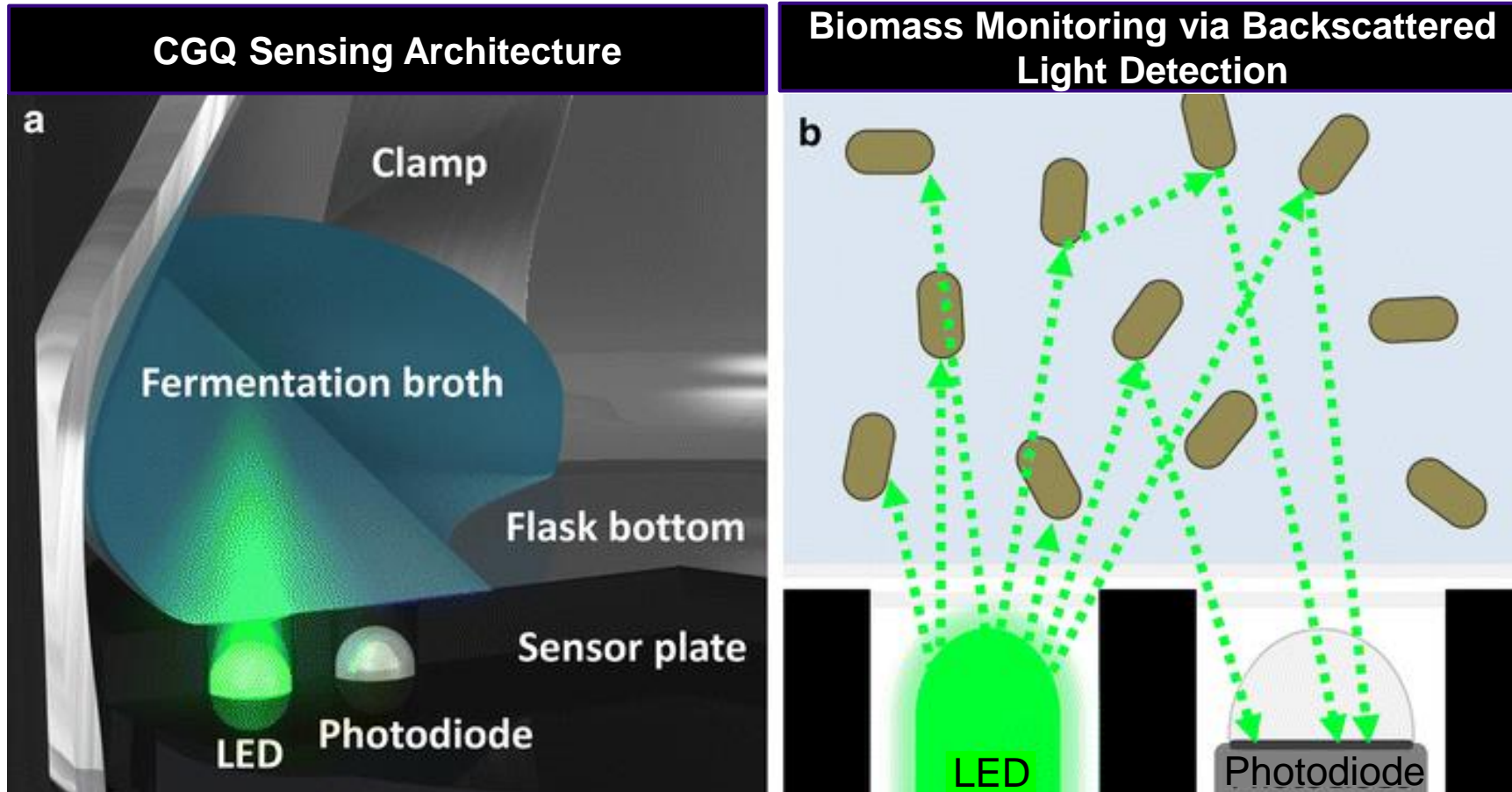


Using different LEDs (green: 521 nm and red: 730 nm), the best wavelength for your phototrophic application can be chosen.

Sensors with Green and Red LEDs

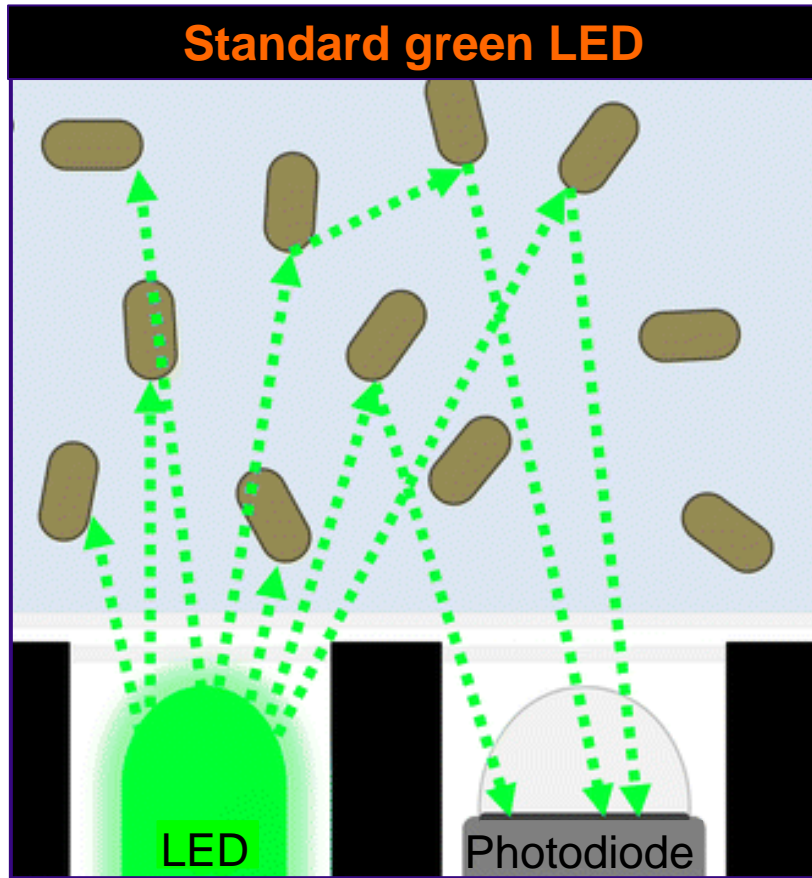


Green LED, 521 nm

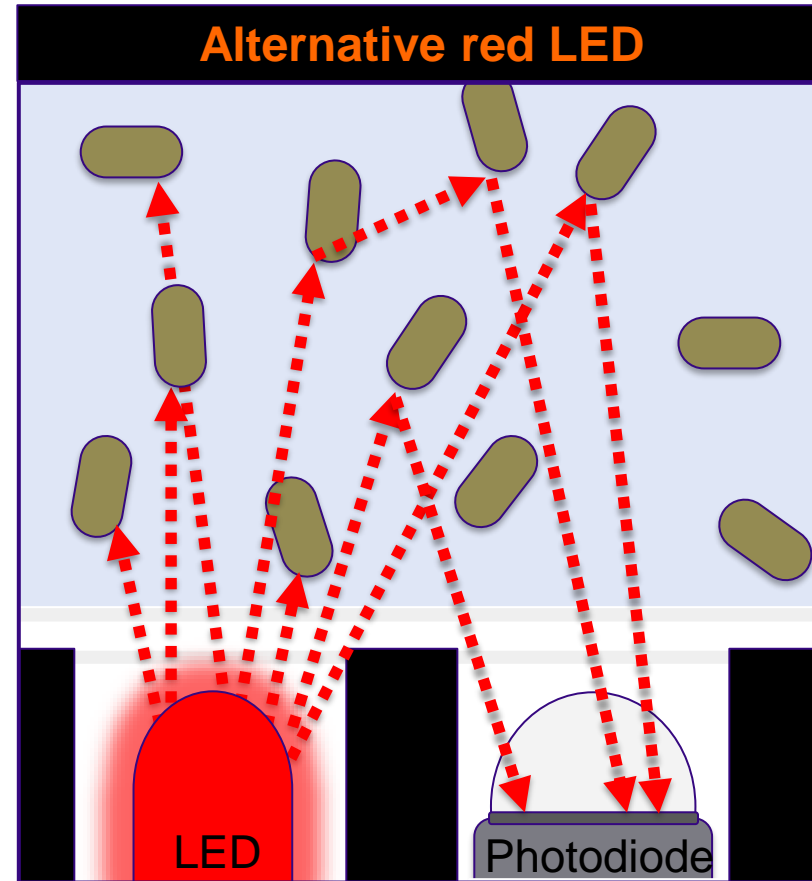


Red LED, 730 nm

For every wavelength, an additional filter can be applied to prevent signal interferences due to too much external light.

Sensors with Red LEDs with and without Filter

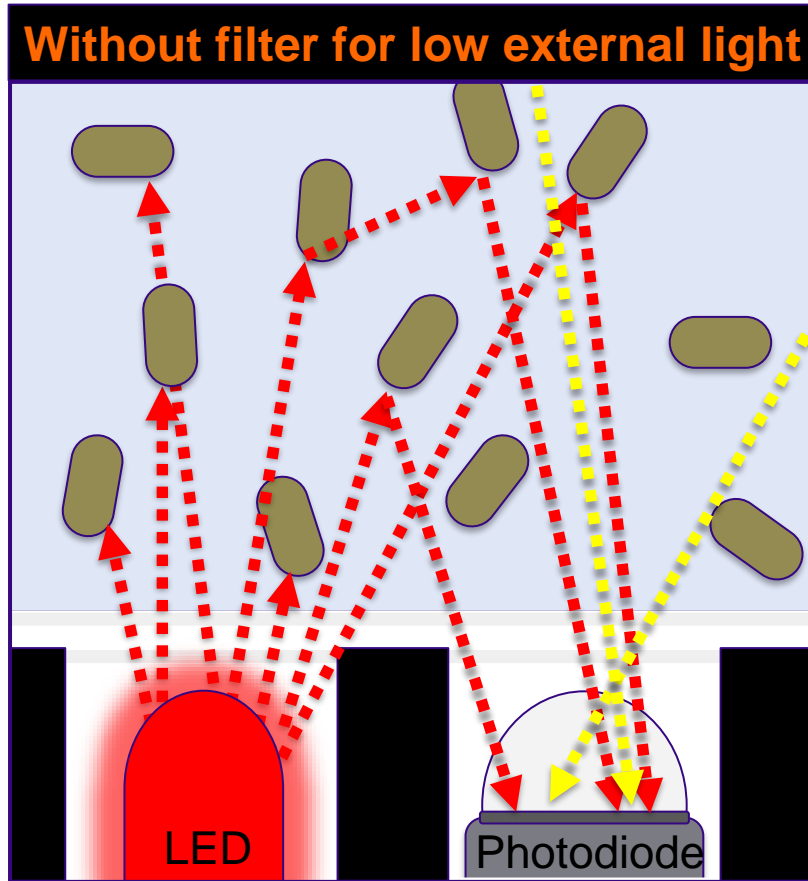





Red LED, 730 nm, no filter

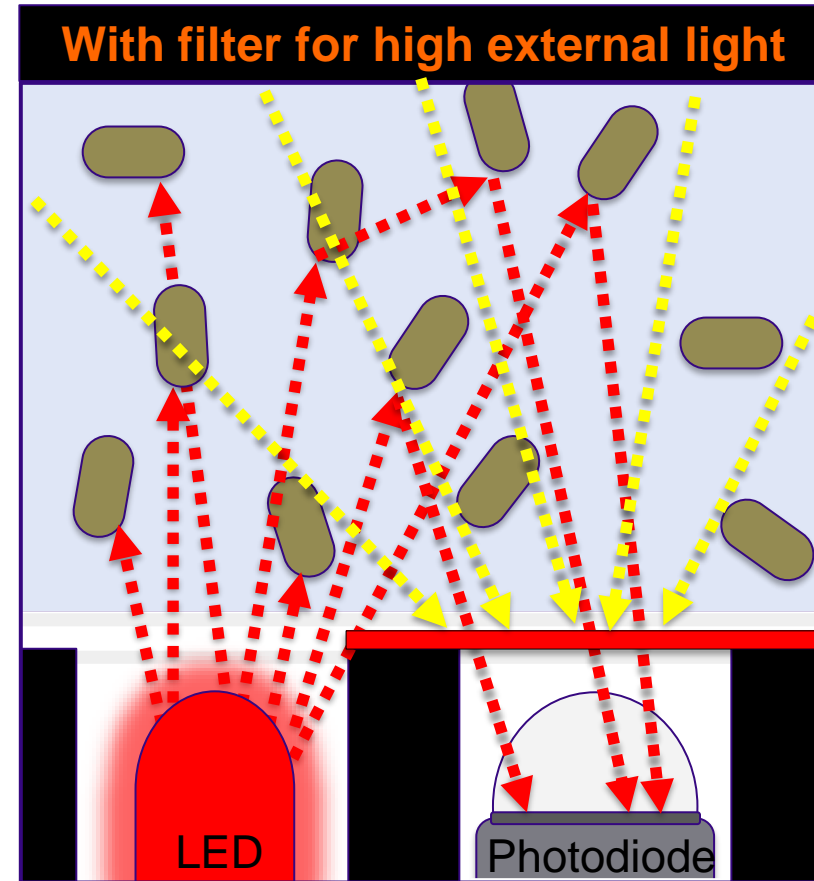

Red LED, 730 nm, additional filter


For both CGQ systems, four different sensors exist for phototrophic trials and can be tested simultaneously.

Different Sensors and Their Specifications

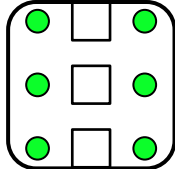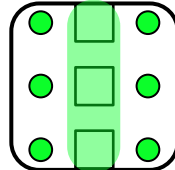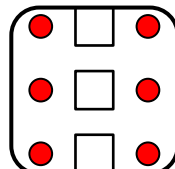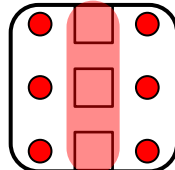LED type	Hardware filter	Schematic	CGQ sensor number	CGQ BioR sensor number
Green LED 521 nm	No Filter		CGQ-SP-F-00198	CGQ-SP-BC-0004
Green LED 521 nm	Filter		CGQ-SP-F-00186	CGQ-SP-BC-0005
Red LED 730 nm	No Filter		CGQ-SP-F-00028	CGQ-SP-BC-0004
Red LED 730 nm	Filter		CGQ-SP-F-00029	CGQ-SP-BC-0006

The CGQ can be used for online monitoring of *Chlorella vulgaris* shake flask cultures and shows good correlation with offline data.


CGQ (850 nm, no filter), OD (750 nm) & Cell Dry Weight (CDW) Measurements: *Chlorella vulgaris*

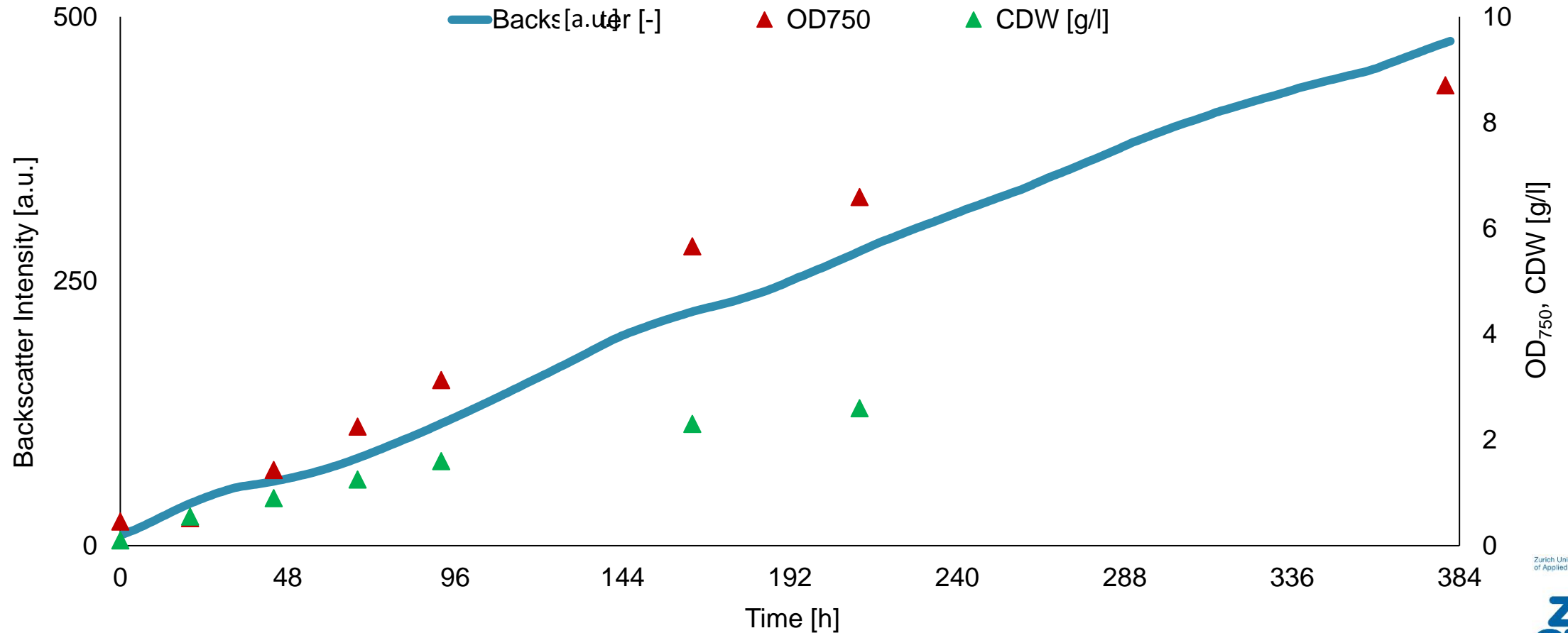

Chlorella vulgaris, 50 mL BBM Medium, 250 mL Shake Flask, 28 °C, 130 rpm

The CGQ BioR is capable of monitoring *Microcystis aeruginosa* in balloon bioreactors and the data correlates nicely with optical density (OD).

CGQ (730 nm) & OD (750 nm) Measurements: *Microcystis aeruginosa* CBT 480

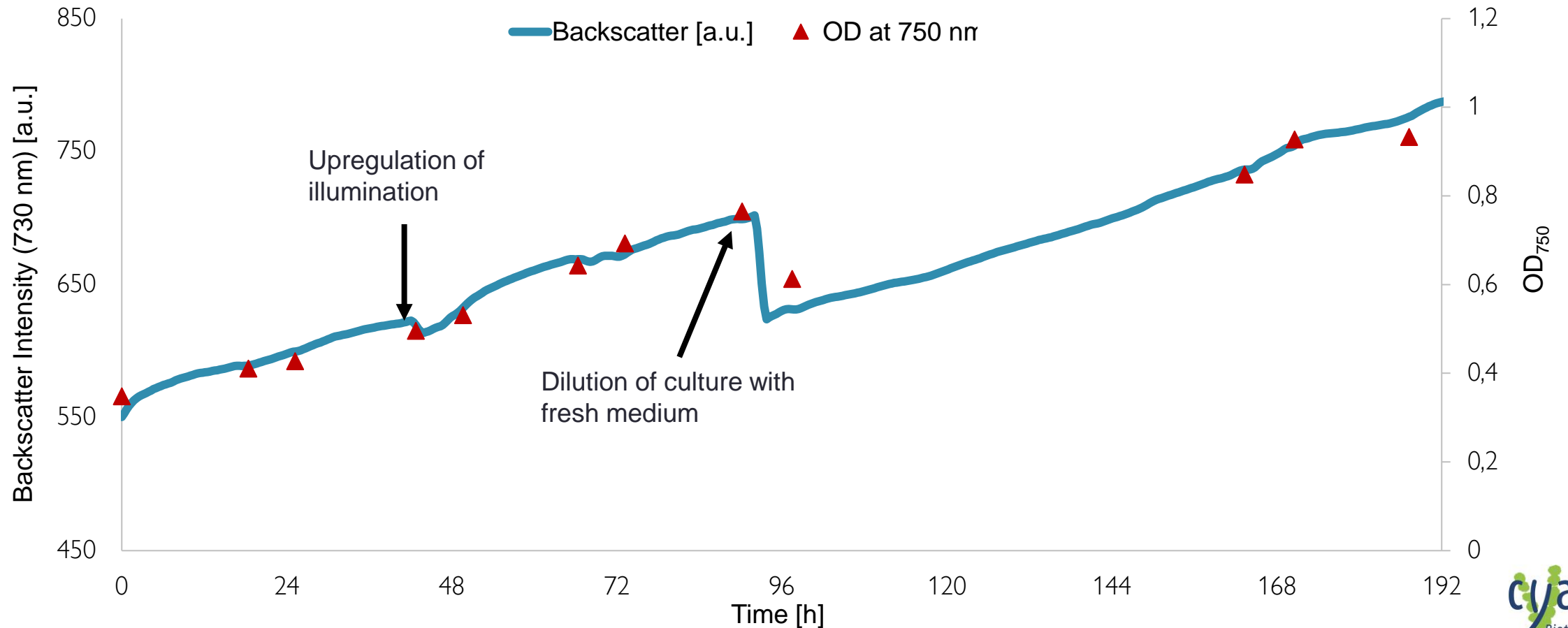


Microcystis aeruginosa CBT 480, 20 L BG11 Medium, 20 L balloon bioreactor, 28 °C



When monitoring *Synechococcus elongates*, the CGQ data correlates nicely with cell dry weight (CDW) and cell number.


CGQ (730 nm+filter), Cell Dry Weight (CDW) & Cell Number Measurements: *S. elongatus*

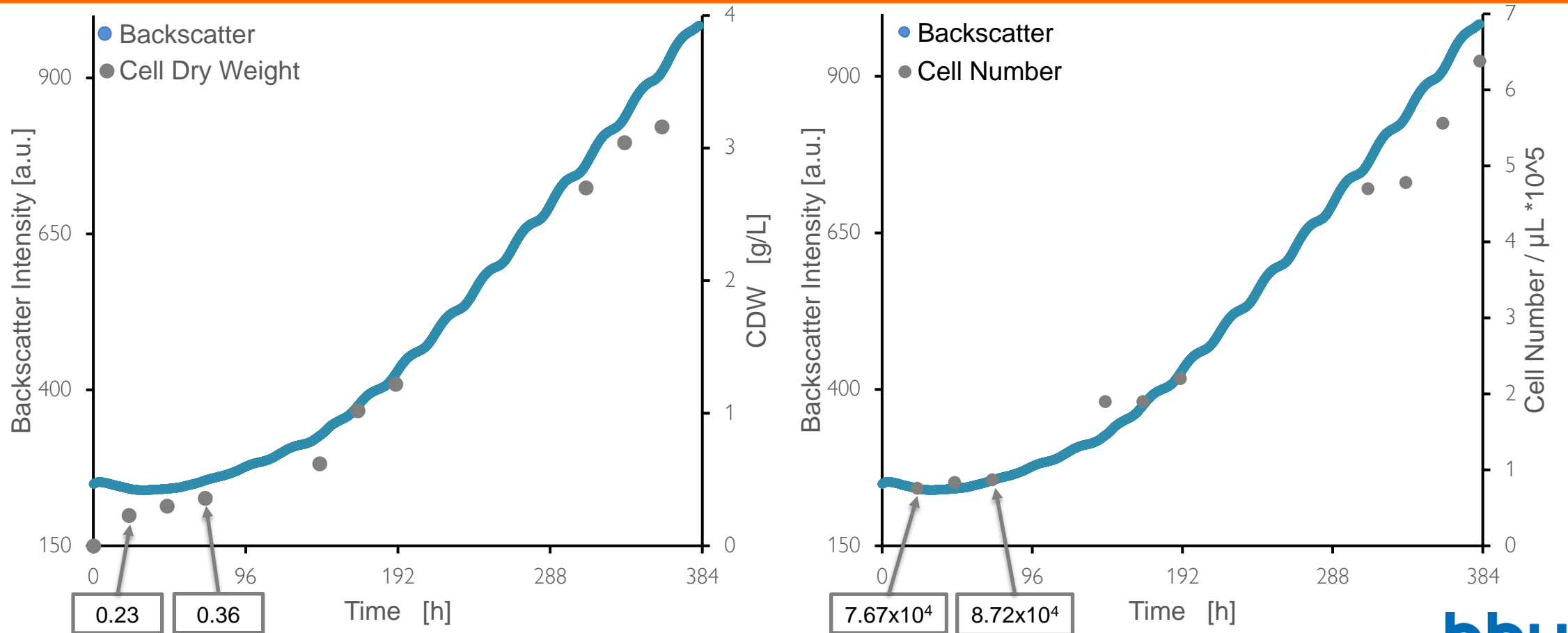

Synechococcus elongates, Medium: BG11, 100 mL in 250 mL flask, 80 $\mu\text{Einstein}$ constant illumination, 150 rpm, 30°C

The CGQ can monitor *Synechocystis sp.* and the data correlates nicely with cell dry weight (CDW) and optical density (OD) at different wavelengths.


CGQ (521 nm, with filter), CDW & OD Measurements: *Synechocystis sp.*

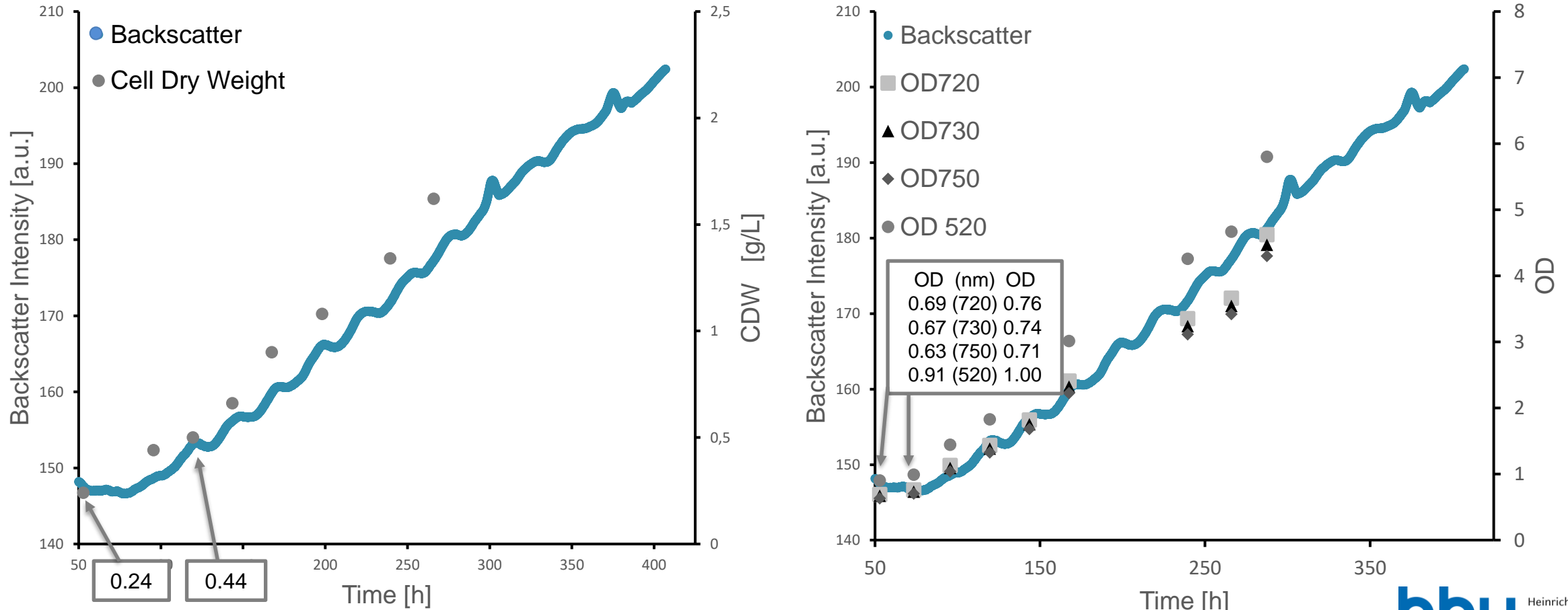

Synechocystis sp., Medium: BG11, 100 mL in 250 mL flask, 80 μ Einstein constant illumination, 150 rpm, 30°C

When monitoring *Arthrospira platensis*, the CGQ data correlates nicely with cell dry weight (CDW) and optical density (OD).


CGQ (521 nm, with/without filter), CDW & OD Measurements: *Arthrospira platensis*

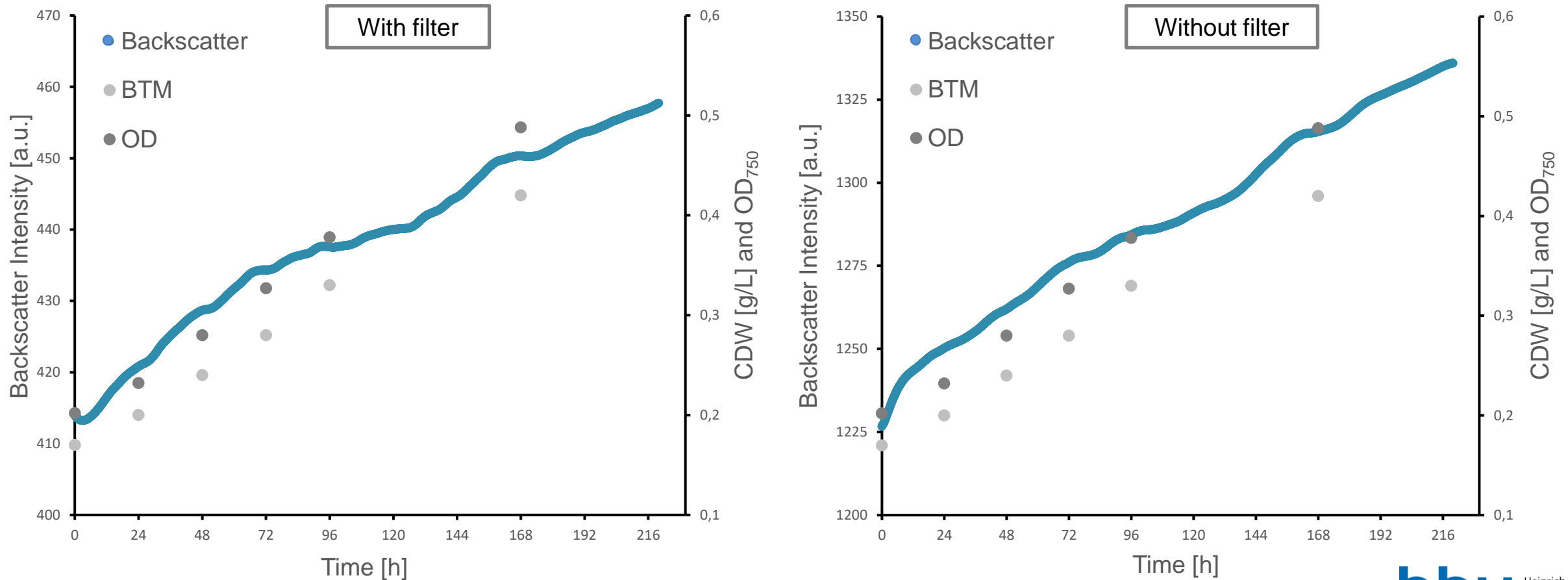

Arthrospira platensis, 100 L Photobioreactor, 150 $\mu\text{mol photons m}^{-2} \text{s}^{-1}$ constant illumination, 2 L/min gassing rate, 26°C

The phototrophic system can be used in a shaker with simulated day-night cycle*.

CGQ Measurements: *Synechocystis sp.* PCC 6803

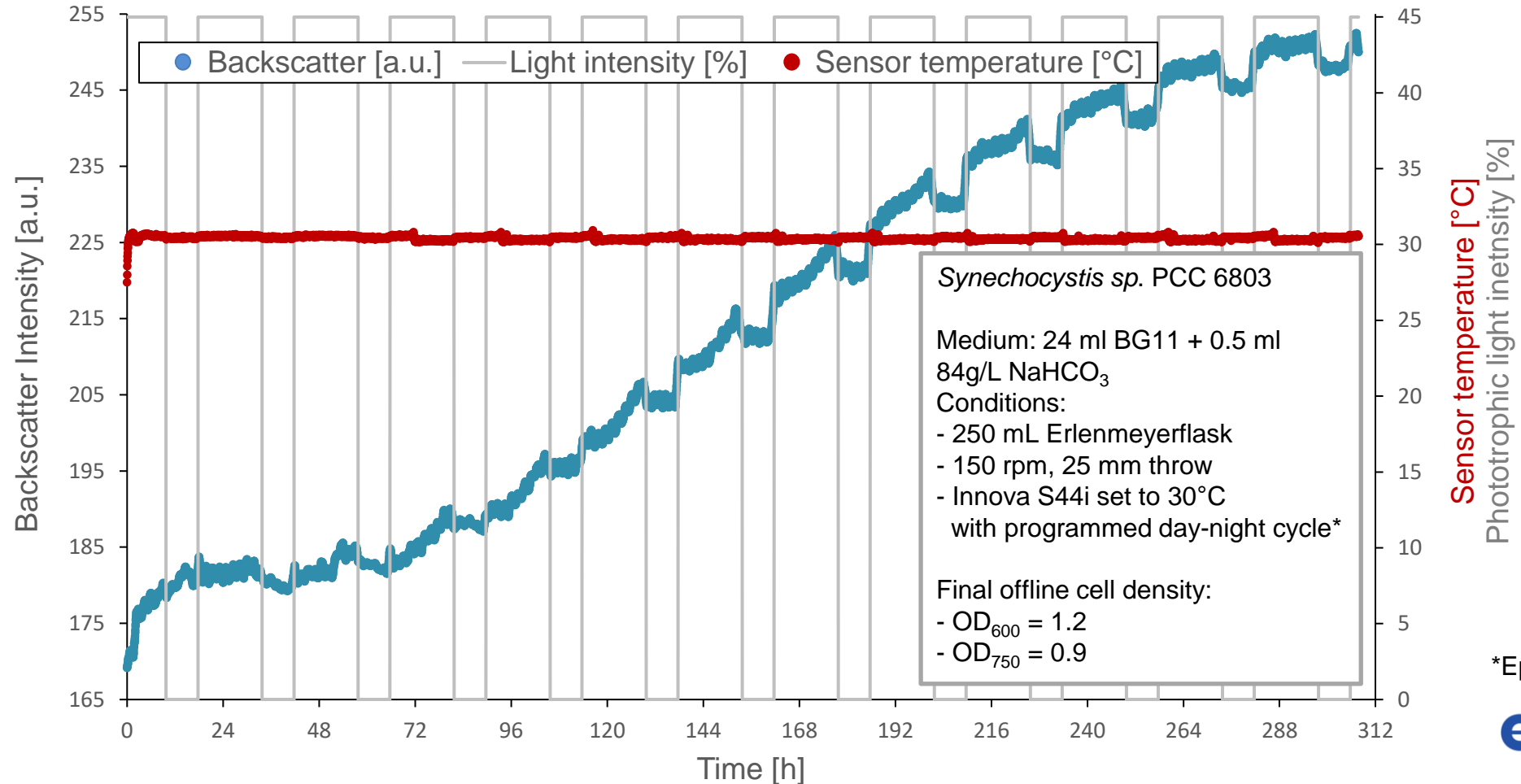

*Eppendorf Application Note

eppendorf

sbi SCIENTIFIC
BIOPROCESSING

Let's Connect!

insights@scientificbio.com
www.scientificbio.com

@ scientific bioprocessing



@scientific bioprocessing



@scientific_bio

